



JELIGHT COMPANY, INC.

2 Mason, Irvine, CA 92618

Phone: (949)380-8774 / Fax: (949)768-9457

E-mail: sales@jelight.com

OZONE GENERATOR MODEL 2001

INSTRUCTION MANUAL

LAMP INSTALLATION

1. Unplug AC power cord from the receptacle on ozone generator.
2. Remove top cover, and lamp access cover from front of unit.
3. Carefully slide lamp assembly into ozone tube, then slide the collar over the lamp wires onto the ozone tube. Make sure the o-ring on the lamp assembly flange is seated around the end of the ozone tube. Secure by tightening.
4. Attach ring terminals from lamp assembly to the power supply. (See page 2, Schematic)

TUBE INSTALLATION

1. The media intake is on the rear of the unit and is marked "MEDIA INLET". Purge oxygen, air, or any another sample you require into this 1/4" tube fitting.
2. The ozone supply is on the front of the unit and is marked "EXHAUST OUTLET" connect this 1/4" tube fitting to its desired destination. Check all fittings for leaks.

Plug in the AC power cord and turn on power switch. Let unit operate for a minimum of five minutes to properly stabilize. Adjust flow gauge to proper setting. The ozone generator is now ready for service.

NOTE:Lamp should be replaced after 5000 hours of use.