



JELIGHT COMPANY, INC.

2 Mason, Irvine, CA 92618

Phone: (949)380-8774 / Fax: (949)768-9457

E-mail: sales@jelight.com

OZONE GENERATOR MODEL 600

I N S T R U C T I O N S

The inlet is located on the side marked "IN". Purge oxygen, air, or whatever sample you require into this end; then purge the sample out through the side marked "OUT".

At the side marked "IN", make sure that the fitting is loosened then slide the sleeve in (open end first) until it stops. There is an O-ring at this spot, therefore you must apply slight pressure to slide it further. Once it completely stops, tighten the fitting snugly.

At the other end, loosen the fitting and slide the lamp into the chamber making sure that the lamp goes into the sleeve. Push the lamp until it completely stops. Once again apply slight pressure to push the lamp through another O-ring; it will completely stop once the lamp is to the end. Tighten the fitting.

At this point, no ozone will be produced. To adjust the ozone concentration, loosen the fitting at the end marked "IN". Slowly slide out the fitting to the desired amount, then tighten the fitting. The ozone rate will change according to your adjustments.